Klean-Contor GmbH

www.lidbag.de
KleanContor@t-online.de

Hotline [49] (0)180 / LIDBAGS
oder [49] (0)180 / 5432247

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The sterilisation bag for highest hygienic requirements – the lid bag for autoclaving

- Easy-to-close, odor- and liquid-tight seal
- Integrated handle makes for easy handling
- Sterilization at temperatures of up to 134°C
- Highly economical and efficient to use
- Meets the most stringent hygienic and occupational-safety requirements
- Suitable for L3 and S3 laboratories as well
- Patented world-wide

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Biologically-contaminated waste is placed in the auto-claving bag right where it is generated. The Desa-Trans container makes the auto-claving bag easy to handle. Opening/closing via foot pedal keeps your hands off the column, thus keeping it uncontaminated. Lid closure is hydraulically cushioned, preventing any unnecessary contamination of the immediate area due to air set in motion (from cover movement).

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Even when the lid of the bag is simply closed (and not sealed), no odors can escape. When the lid bag is full, simply press down the lid to seal it. Once sealed, the bag is odor-tight and leak-proof. Due to the patented sealing process, no odors can escape during the actual sealing of the bag.

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After the bag is sealed, it can be easily removed from the Desa-Trans column using its integrated handle and replaced by a new lid bag. The full and sealed bag will be taken off to be autoclaved on the premises. In most cases, the lid bag will not require any additional container for transporting it.

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Using the integrated handle the lid bag is placed in the autoclave. The lid bag and its contents are then disinfected or sterilized at either 121°C or at 134°C.

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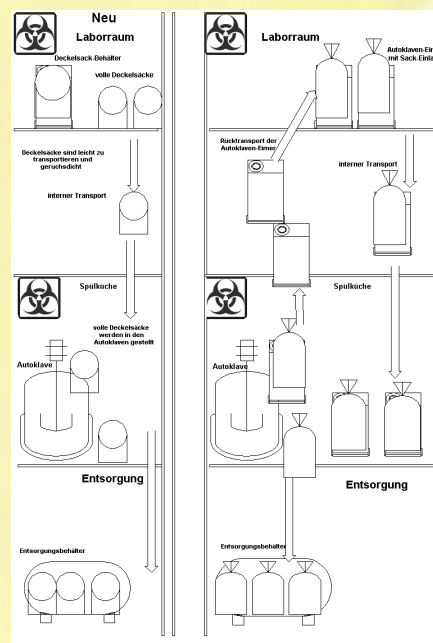
After the autoclave has cooled, the lid bag can be taken out. Any liquid components cannot escape from the lid bag after the autoclaving either. This means that transporting the lid bag to the disposal site is clean and hygienic. The key logistical advantage: because additional containers are not necessary for transporting the lid bag on your premises, there are no additional containers to be disinfected and transported back to where the waste is generated.

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The lid bag – a highly-economical and efficient solution for dealing with bio-hazardous waste. The lid bag also meets the highest hygiene and labor-safety regulations – even for L3 and S3 laboratories.

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- odor- and liquid-tight
- adhesive seal
- liquid binding agent
- for high safety standards

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- BAM certification
- sturdy and puncture-proof
- odor- and liquid-tight
- integrated handle

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- high-grade, polished stainless steel
- comes with non-skid circular stand or with 'roll boy' (rollers)
- available for lid bags with 6 - 60 liter capacity

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- hygienic handling
- hygienic storage
- odor-tight
- leak-proof
- no refrigeration required